



## NEW features dataLink / dataLink Pro !

### Network connection


Network connectivity has been improved. Just like a standard printer, the dataLink **can now be connected to a network**. This allows printing on one or more cab printers **without needing a USB cable**, and from one or multiple computers. In this way, **print configurations are memorized even when switching users**.

### Label Information and Formatting

A **warning message** now notifies the user if the code is too long to fit on the label. The maximum length for the code to fit is **73 alphanumeric characters**. This limit may be reduced if the code includes **special characters** or **uppercase letters**. This rule applies to both the **easySpiral Management** module and the **Label Printing** module.

### Automatic Import

It is now possible to configure an **automatic import of .csv files**. In practice, the user saves a template Excel file in a defined location, which is then linked in the software. The operator updates the file at this usual location before each print session, and the import into the printing modules (**easySpiral Management** and **Label Printing**) happens automatically each time the module is opened.

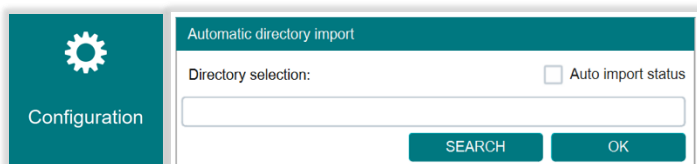

### Print Options

- The system now supports **QR code** printing.



- Print size adapts automatically**: if the content fits on a **single line**, the font size is increased.




- A **"symbol" column** has been added. To add a specific symbol, the user refers to a table where codes are associated with the available symbols. Users can customize symbol codes in the **Configuration module** (ex: ■ = Total flora, ▲ = yeast / mold).
- Rows of printed labels are now **highlighted in blue**.
- From now on, **matte labels** will be used.

Symbol	Value
●	total_cnt
▲	yeast_mold
◆	e.coli
■	4
▨	5
□	6

### User Management

- LDAP** management is now available.
- Automatic saving** is no longer managed in the account management module. It is now accessible in the **Configuration module**. This makes it possible to set up automatic saving even if account management is not activated.

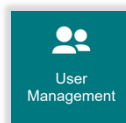

☐ Local configuration ☒ LDAP configuration

☐ Ldaps

URL

Root domain (base DN)

Login

Password

Timeout (in ms)