

# BagFilter Pull-Up®

*Lateral filter bag with pull up system*

The BagFilter Pull-Up® has been especially designed to make your pipeting easier with any type of pipet and very short tips ( $\geq 4$  cm). With a unique and patented filter Pull-Up® system, just pull up the filtrate for easy pipeting. The BagFilter Pull-Up® is a “must” for accurate sample preparation and special tests such as PCR.


**Accurate:** ideal for very precise pipeting ( $\leq 5 \mu\text{L}$ )

**Adaptable:** can be used with very short tips

**Economical:** same bag is used to blend, filter and pipet

**Sterile:** eliminates cross contamination

**Practical:** can be closed with a BagClip® or heat-sealed

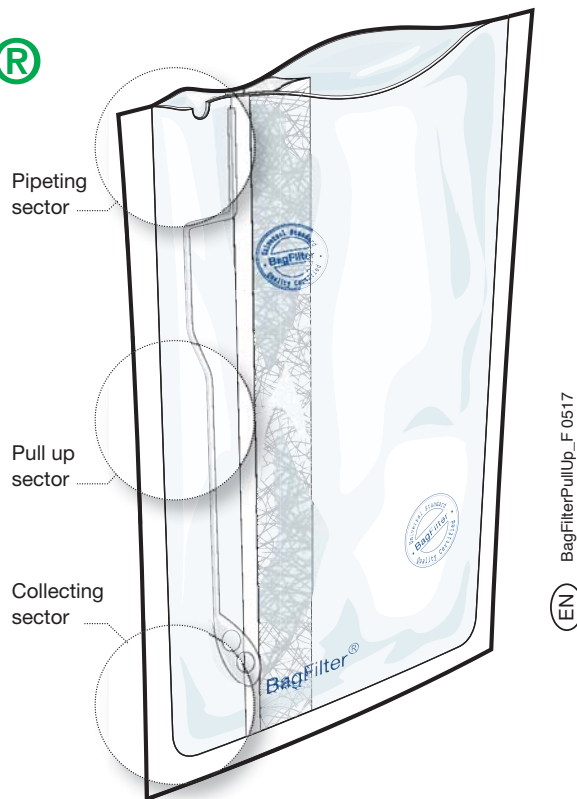

[www.interscience.com](http://www.interscience.com)

**interscience INTERNATIONAL**

30, ch. Bois Arpents - 78860 St Nom - FRANCE  
T: +33 (0)1 34 62 62 61 - F: +33 (0)1 34 62 43 03  
Email: [info@interscience.com](mailto:info@interscience.com)

**interscience USA & CANADA**

32 Cummings Park - Woburn, MA 01801 - USA  
T: +1 781 937 0007 - F: +1 781 937 0017  
Email: [sales.usa@interscience.com](mailto:sales.usa@interscience.com)

**interscience CHINA**

上海徐汇区吴兴路277号锦都大厦588室-200030  
电话: +86(0)21-64739390 - 手机: +86 189 3097 0733  
邮址: [sales.china@interscience.com](mailto:sales.china@interscience.com)


**interscience SOUTH-EAST ASIA**

541 Orchard Rd - 09-01 Liat Towers - SINGAPORE 238881  
T: +65 6933 1389 - M: +65 8549 1217  
Email: [sales.asia@interscience.com](mailto:sales.asia@interscience.com)

# Fast, accurate and easy


## How to use the BagFilter Pull-Up®

By pulling up the liquid up, it becomes simple to pipet a small part of the filtrate. **1** - Just pinch the bag between thumb and index fingers to catch the filtrate into the pull up sector. **2** - Transport the liquid up. The filtrate is temporarily fixed in the pipeting pocket. **3** - You can now pipet with very short tips ( $\geq 4\text{cm}$ ) and after a while, the pocket empties itself by gravity.

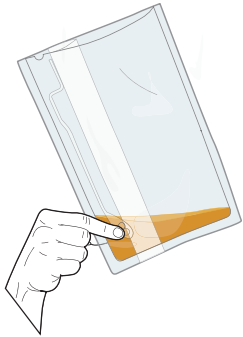


### 1 - Pinch

Pinch the bag to pull up the liquid.

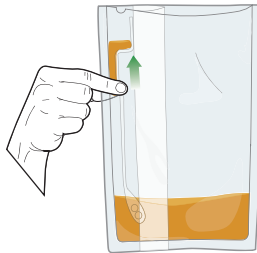

### 2 - Pull up

Pull up the filtrate to the top of the bag.

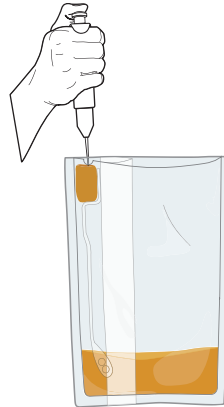

### 3 - Pipet

Pipet easily the filtrate in the pipeting pocket.

### Pull-Up® PATENTED

Filter pump system created by interscience for easy pipeting with short tips.



Pull up sector

## Technical specifications

- **Size** : 190 x 300 mm
- **Capacity** : 50 - 400 ml
- **Pocket volume** :  $\approx 1000 \mu\text{l}$
- **Multilayer® film**
- **Lateral non woven filter** : 50  $\mu\text{m}$
- **Minimum tips length** : 4 cm
- **Quantity per box** : 500

Ref : 111 625