

GeckoGrip for BagOpen®

Bag attachment system (Patented) on DiluFlow®
Ref. 505 009 (400/3500 mL) - **Ref. 505 015** (80/100 mL)



Manufactured:



This patented system replaces adhesive rubber and keeps the test bag glue-free, even on wet surfaces. The design of GeckoGrip strips guarantees splash-proof and detergent-proof adhesion, as well as integrity that eliminates the risk of cross-contamination.

OPERATING MODE

1. Center and stick a GeckoGrip strip on the inside of each end of the BagOpen®.
2. Slide a new bag between the BagOpen® walls.
3. Let the bottom of the bag rest on the base of the accessory.
4. Pinch the moving ends of the BagOpen® until the GeckoGrip strips fully adhere to the sides of the bag.
5. Release to open the bag.

MAINTENANCE

1. Apply a little soap to hands. Soap the surface of the GeckoGrip strips.
2. Lather under cold water. Rinse off.
3. Dry BagOpen® walls with paper towels.
4. Dab the surface of each GeckoGrip strip with the paper towel.
5. Pour 70% ethanol over each GeckoGrip strip to disinfect.
6. Remove excess alcohol by dabbing each GeckoGrip strip with the paper towel.
7. Replace BagOpen® on dilutor. Place the bag according to operating mode.

TECHNICAL SPECIFICATIONS

Different GeckoGrip models are available, depending on bag dimensions and the type of BagOpen® mounted on DiluFlow®.



GeckoGrip Ref. for BagOpen®	Ref. 505 009	Ref. 505 015
Compatibility with BagOpen®	Ref. 211 040 (400 mL - 2000mL) Ref. 211 350 (3500 mL)	Ref. 211 010 (80/100 mL)
Material	Silicone rubber	
Manufacturing process	Injected film	
Number of surface adhesion points	29000 /cm ² approx.	
Weight	300 g/m ² approx.	
Color	natural	
Storage conditions	-50 °C - +150 °C approx. / -122 °F - +302 °F approx.	
Packaging	Pack of 10 adhesives	
Belt thickness	Approx. 0.35 mm	
Bandwidth	25 mm	
Belt length	90 mm	40 mm
Finishes	Rounded corners	Cut edges