

**NEW**

# *insta***BAG**<sup>®</sup> MILK

Filter blender bag with dehydrated media



Your chocolate analyses improved!

*insta***BAG**<sup>®</sup> MILK is a ready-to-use sample preparation bag for instant chocolate product analysis.

**interscience**

# Ready-to-use media for easy preparation



## Just add warm water

The pre-dosed milk powder bag is directly integrated into the bag. Simply add a sample and warm purified sterile water. The milk powder dissolves completely during mixing. Preparation is immediate.

The pre-dosed dehydrated medium is French milk powder from Normandy. *instaBAG*<sup>®</sup> MILK can therefore be used for the analysis of chocolate products: cocoa powders, chocolate bars and various finished chocolate products.

1

Place the sample in an *instaBAG*<sup>®</sup>



2

Add sterile and warm deionized water



3

Blend the *instaBAG*<sup>®</sup> with the BagMixer<sup>®</sup>



**The sample is ready to be pipetted**

*instaBAG*<sup>®</sup> MILK is available in **225**, **900**, **2250** and **3375 g**.

**Request free samples!**

Contact us for **Infos | Videos | Demos | Quote** at [info@interscience.com](mailto:info@interscience.com)

and visit [www.interscience.com](http://www.interscience.com)